LabSpeed – Table Export to SQL Server Database User Guide

Overview

The **Table Export SQL Server extension** adds the ability to export data from a LabSpeed Table View directly to a SQL Server database.

The LabSpeed Table View provides a drag & drop mechanism for creating a flat table from relational instrument data. The columns in this table represent any subset of the available instrument data and may be arranged in any order. In addition, the table may be quickly rearranged in any of three distinct styles. Once setup, the data may be exported as Delimited Text, Excel (.xls) or to a SQL Server database table. Exporting to a SQL Server database requires an extension add-in be installed.

Add-in Installation

The Table Export to SQL Server LabSpeed add-in installer may be obtained by visiting the Topos Technologies website at <u>www.topostech.com</u> and clicking on the <u>Feature</u> <u>Extensions</u> link. Download and run the <u>ToposTableExportSQLServer.msi</u> installer program and follow the directions.

When LabSpeed is next run, you will be directed to the Client Manager in order to add the new Table Export SQL Server Feature Add-In to LabSpeed's client list.

SQL Server DB Connection

The SQL Server connection parameters must be setup correctly. Ensure parameters are correct by choosing the menu *Tools->Client Connections*. This is where all the connections for all Clients are managed.


Select the Client "Topos : TableExportSQLServer" and the connection named "Export Table to SQL Server" and click *Edit*. Or create a new connection. Enter all required SQL Server log on fields, including Server Name, User Name and Password and then select a Database. Test the connection to makes sure it is valid and then save the connection to the same name (which overwrites the existing connection) or to a different name. If you save the connection to a different name, remember to make it the *Active Connection*.

Table Export SQL Server Setup

While in Table Designer view, click the Export toolbar button. The Table Designer Export dialog is where you choose the SQL Server Export provider and setup column maps.

The Table Export SQL Server Provider

On the General tab, select the Export to SQL Server export provider. If it does not appear in the drop down list, it may need to be installed. The Export to SQL Server option is a special LabSpeed Add-In feature.

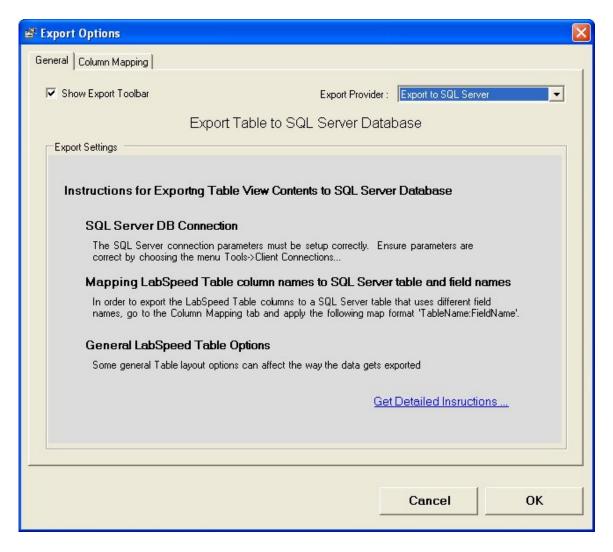

General instruction appear describing the process of setting up for exporting to SQL Server. In addition, a button is available for obtaining more detailed instructions.

Mapping Table column names to SQL Server field names

Each LabSpeed Table column that you wish to export must be mapped to the corresponding SQL Server table column name.

On the Column Mapping tab, enter field names that map the Table column name to the SQL Server table column name by applying the following rules:

- 1. Map format TableName:FieldName. Leading and trailing spaces are not important and will be removed
- 2. Any mapped items that are not of the TableName:FieldName format will be ignored.
- 3. The same Table Name has to be used for all mapped items

For example, if the SQL Server database table name is "Sample", the following maps might be applied:

Export Options				
General	Co	olumn Mapping		1
8	<u> </u>	Field Name	Map Name	
	1	SampleName	Sample:SampleName	
		AcquireDate	Sample:Date	
		SampleType	Sample:SampleType	
		User Name	Sample:Operator	
	1	SampleWeight		
		InitiaMolume		
		FinaVolume		-
		CorrectionFactor	Sample:Factor	1
		As1890	Sample:As	
		Ca3179	Sample:Ca	
		Ca3933		1
		Cu3247	Sample:Cu	
		Hg1849	Sample:Hg	
		K_7664	Sample:K	1
		Mg2852	Sample:Mg	
		Na5889	Sample:Na	1
		Na5895		1
		Рь2203	Sample:Pb	
	Γ	F Show All Maps	Field Type:	3
			Cancel	ок

Other Mapping Rules

- 1. The table and mapped fields must exist in the SQL Server Database. It does not matter if there are additional, unmapped database fields in the SQL Server Table.
- 2. The Database column types must match the associated LabSpeed field data type. For example, if the LabSpeed data type is "Int32", then the SQL Server table field must also be "Int32". However, any LabSpeed field type may be exported to a database string field. Click on the Field Types... button to view the LabSpeed exported fields and their Field Types.
- 3. If you wish to turn off the Table Option "Show duplicate primary labels", the primary columns in the Database must have the "allow null" column option checked.

Note: Changing the Table Style (*Standard*, *Value Type* and *Value Type Column*) will change the format of the table and result in more or fewer columns exported. This will affect maps as well.